



# SONOCHECK® ABD06

Non-contact Bubble Detectors | Bioprocessing

SONOTEC 

Process monitoring of liquids in biopharma lab and production environments enables continuous control and optimization of critical process parameters. Hence, the early detection of air bubbles is an effective instrument for making processes more stable, productive and efficient, and reducing costs in the long term. Highly precise and flexible **non-contact clamp-on SONOCHECK® ABD06 bubble sensors** are perfect for continuous air bubble monitoring and full/empty detection.

SONOTEC® implemented a patented closed loop algorithm to ensure a fast detection of even smallest bubbles and to guarantee constant bubble sensitivity widely independent of the quality of the acoustic coupling. SONOCHECK® ABD06 bubble sensors can spot bubbles down to 1/3 of the inner diameter of the tubing. Thus, the bubble detectors profit from SONOTEC's long-term experience in developing medical grade bubble detectors for hemodialysis and ECMO machines.

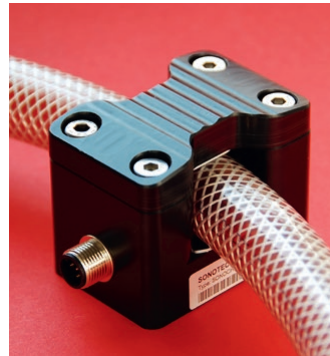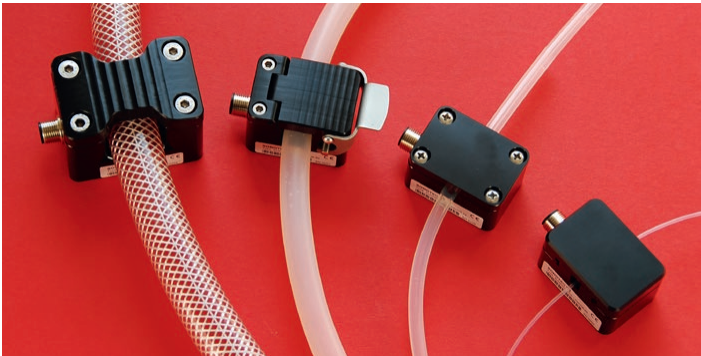

## KEY FEATURES

- High grade sensors for air bubble monitoring and full/empty detection
- Compact sensor design with integrated electronics for data processing
- Non-contact clamp-on mechanism enabling direct measurement through the tubing wall
- For a wide range of bioprocessing-specific plastic tubing, such as platinum cured silicone, TPE, reinforced tubing, PVC, PE, PFA, FEP
- No contamination and no shear stress on cells due to contact-free sensor
- Suitable for WFI, cell culture media, buffer solutions, and most aqueous solutions
- Stable sensor reliability to changing acoustic conditions, such as moderate changes in buffer solution concentration, fluid density as well as viscosity
- Optional sensor version with ATEX certification for use in hazardous areas

## APPLICATIONS

|                                   |   |
|-----------------------------------|---|
| <b>BIOREACTORS / FERMENTATORS</b> | Detection of excessive foaming in feed/harvest lines  |
| <b>CHROMATOGRAPHY</b>             | Prevention of air being pumped into columns & diversion of air infused liquids around the columns |
| <b>TANGENTIAL FLOW FILTRATION</b> | Prevention of air entry into filter cassettes   |
| <b>FEED STREAM</b>                | Continuous monitoring for air bubbles in the feed to prevent from air entry into the filter       |
| <b>PUMP PROTECTION</b>            | Detection of air bubbles caused by cavitation and protection from dry running                     |
| <b>TRANSFER LINES</b>             | Notification, when transfer reservoirs run dry  |
| <b>FILL &amp; FINISH</b>          | Detection of undesired air entering vessels to be filled  |



# SONOCHECK® ABD06 | TUBING SIZE CHART

| Tube OD<br>inch | Soft / Flexible Tubing |                  | Hard Tubing<br>PE, PTFE, FEP, PFA, Teflon |                  | Braided Tubing       |                  | RS485 | UART | 4-20 mA | PNP |
|-----------------|------------------------|------------------|---|------------------|----------------------|------------------|-------|------|---------|-----|
|                 | SONOCHECK®<br>Sensor   | Lid<br>Order No. | SONOCHECK®<br>Sensor                      | Lid<br>Order No. | SONOCHECK®<br>Sensor | Lid<br>Order No. |       |      |         |     |
| 0.0625          |                        |                  | ABD06.85                                  | Integrated       |                      |                  |       | ✓    |         | ✓   |
| 0.125           |                        |                  | ABD06.100                                 | Integrated       |                      |                  |       | ✓    |         | ✓   |
| 0.250           | ABD06.115*             | 200 03 0009      | ABD06.115*                                | 200 03 0010      |                      |                  | ✓     |      | ✓       | ✓   |
| 0.375           | ABD06.120              | 200 03 0055      | ABD06.114*                                | 200 03 0035      |                      |                  | ✓     |      | ✓       | ✓   |
| 0.438           | ABD06.120              | 200 03 0060      |   |                  |                      |                  | ✓     |      | ✓       | ✓   |
| 0.500           | ABD06.117              | 200 03 0054      | ABD06.117                                 | 200 03 0049      | ABD06.117            | 200 03 0054      | ✓     |      | ✓       | ✓   |
| 0.563           | ABD06.117              | 200 03 0052      |   |                  |                      |                  | ✓     |      | ✓       | ✓   |
| 0.625           | ABD06.117              | 200 03 0060      | ABD06.121                                 | 200 03 0054      | ABD06.117            | 200 03 0057      | ✓     |      | ✓       | ✓   |
| 0.750           | ABD06.121              | 200 03 0058      | ABD06.121                                 | 200 03 0057      |                      |                  | ✓     |      | ✓       | ✓   |
| 0.875           | ABD06.121              | 200 03 0061      |   |                  | ABD06.121            | 200 03 0061      | ✓     |      | ✓       | ✓   |
| 1.000           | ABD06.102              | Integrated       | ABD06.102                                 | Integrated       | ABD06.102            | Integrated       | ✓     |      | ✓       | ✓   |
| 1.125           | ABD06.116              | Integrated       |   |                  | ABD06.116            | Integrated       | ✓     |      | ✓       | ✓   |
| 1.375           | ABD06.123              | Integrated       |   |                  |                      |                  | ✓     |      | ✓       | ✓   |
| 1.405           |                        |                  |   |                  | ABD06.123            | Integrated       | ✓     |      | ✓       | ✓   |


\* Sensors have a screw lid

## MEASUREMENT PRINCIPLE

✓ **SONOCHECK® ABD06 non-contact clamp-on bubble sensors** incorporate intelligent ultrasonic transmission technology. The presence of air bubbles and obstructions is detected by means of dynamic amplitude monitoring.

Ultrasound waves are mechanical waves and are thus subject to the laws of their physics. Depending on the sound impedance of the adjacent media, reflection and transmission take place at the interfaces. If the impedance differences of the adjacent media are small, a transition takes place. At larger differences the sound wave is reflected and does not penetrate the adjacent media.

When an air bubble passes the sensor channel, the signal level drops. Hence, the higher the drop of the signal level, the larger the bubble size.

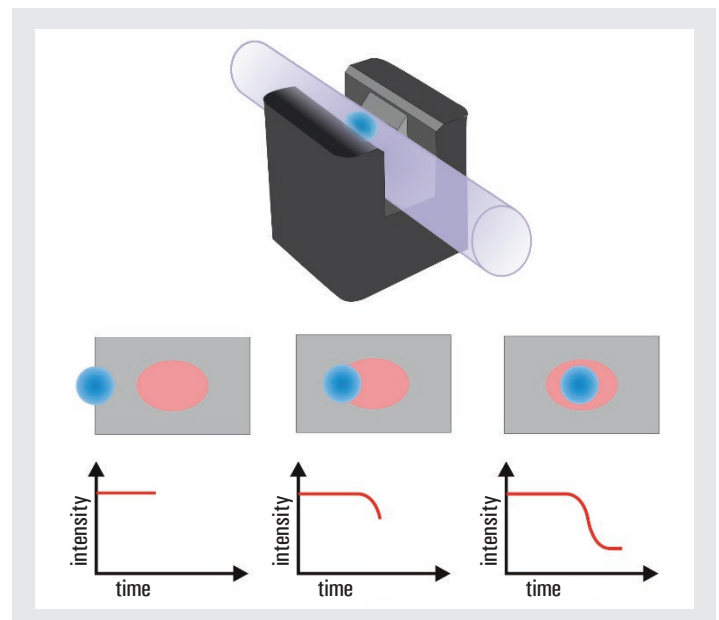

SONOTEC® preserves the right to change technical specifications without further notice. (Rev. 1.1 / 2020-06-19)

### SALES & SUPPORT

**SONOTEC GmbH**  
 Nauendorfer Str. 2  
 06112 Halle (Saale)  
 Germany

**phone** +49 (0)345 / 133 17-0  
**e-mail** sonotec@sonotec.de  
**web** www.sonotec.eu

**SONOTEC** 

Certified according to ISO 9001 & EN ISO 13485